

Division of Signal Transduction Therapy

Standard Operating Procedure

Preparation of active Phosphatidylinositol-5-phosphate 4-kinase type II alpha

Enzyme description:- PIP5K2A [1 - 406]

Clone number:- DU 12296

Source:- Recombinant

Expression system:- *E.coli*

Tag:- N-terminal His(6)

Purification method:- Ni²⁺-NTA agarose

Expression level:- 10 mg/L

Calculated molecular mass:-

Monoisotopic 49,737.04 daltons

Average Mass 49,768.50 daltons

[cysteines reduced, methionines have not been oxidised]

Theoretical pI:- 6.93

Purity:- >80 %

Activation protocol:- Constitutively active

Enzyme storage buffer:-

50 mM Hepes/NaOH pH7, 150 mM NaCl, 5 mM DTT, 20% Glycerol,
1 mM Benzamidine, 0.2 mM PMSF

Storage temperature:- -70 °C [Long term stability to be determined]

Assay:- Kinase Glo

Assay buffer:-

50 mM Tris, 134 mM KCl, 2 mM DTT, 1 mM EGTA, 10 mM MgCl₂,

Substrate:-

PI(5)P diC8 Final concentration: 50 μM

Division of Signal Transduction Therapy

Clone Data Sheet

PIP5K2A [1 - 406]

<u>Protein</u>	PIP5K2A [1 – 406]
<u>Clone number</u>	DU 12296
<u>Species</u>	Human
<u>Accession number</u>	NM_005028.4
<u>Tags</u>	N-terminal His(6)
<u>Bacterial expressed protein</u>	MGSSHHHHHSSGLVPRGSHMASMTGGQQMGRGSMATPGNLGSSVLASK TKTKKHFVAQKVKLFRASDPLLSVLMGVNHSINELSHVQIPVMLMPD DFKAYS KIKVDNHLFNKENMP SHFKFKEYCPMVFRNL RERFGIDDQDFQ NSL TRS APLPNDSQARSGARFH TSYDKRYI IKTITSEDVAEMHNILKKY HQYIVECHGITLLPQFLGMYRLNV DGVEIYVIVTRNVFSHRLSVYRKYD LKGSTVAREASDKEAKELPTLKDNDFINEGQKIYIDDDNNKKVFLEKLK KDVEFLAQQLKLM DY SLLVG IHDVERAEQEEVECEENDGE EEGESDGTHP VGT PPDSPGNTLNSSPPLAPGEFD PNIDVYG IKCHEN SPRKEV YFMAII DILTHYDAKKAAHA AKTVKH GAGAEISTVNPEQYSKRFLDFIGHILT
<u>Native sequence</u>	Amino acids M1 – T406 (end) of human PIP5K2A. Residue M35 of the fusion protein is equivalent to M1 of the native enzyme. The His(6) tag is located at residues 5 – 10.
<u>Protease cleavage</u>	Thrombin (<u>LVPRGS</u>) at residues 14 - 19
<u>Cloning sites</u>	<i>Bam</i> H1 and <i>Not</i> I sites of pET28a

Division of Signal Transduction Therapy

Complete
nucleotide
Sequence

ggatccATGGCGACCCCCGGCAACCTAGGGTCCTCTGTCTGGCGAGCA
AGACCAAGACCAAGAAGAAGCACTTCGTAGCGCAGAAAGTGAAGCTGTT
TCGGGCCAGCGACCCGCTGCTCAGCGTCCTCATGTGGGGGTAAACCAC
TCGATCAATGAACTGAGCCATGTTCAAATCCCTGTTATGTTGATGCCAG
ATGACTTCAAAGCCTATTCAAAAATAAAGGTGGACAATCACCTTTTAA
CAAAGAAAACATGCCGAGCCATTCAAGTTAAGGAATACTGCCGATG
GTCTCCGTAACCTCGGGAGAGGTTGGAATTGATGATCAAGATTTCC
AGAATTCCCTGACCAGGAGCGCACCCCTCCCCAACGACTCCCAGGCCCG
CAGTGGAGCTCGTTTCACACTTCCTACGACAAAAGATACTCATCAAG
ACTATTACCACTGAAGACGTGGCCAAATGCACAAACATCCTGAAGAAAT
ACCACCACTACATAGTGGAAATGTCATGGGATCACCCCTCTCCCCAGTT
CTTGGGCATGTACCGGCTTAATGTTGATGGAGTTGAAATATATGTGATA
GTTACAAGAAATGTATTCAAGCCACCGTTGTCTGTATAGGAAATACG
ACTTAAAGGGCTCTACAGTGGCTAGAGAAGCTAGTGACAAAGAAAAGGC
CAAAGAACTGCCAACTCTGAAAGATAATGATTCAATTATGAGGGCCAA
AAGATTATATTGATGACAACAACAAGAAGGTCTTCCTGGAAAAACTAA
AAAAGGATGTTGAGTTCTGGCCAGCTGAAGCTCATGGACTACAGTCT
GCTGGTGGATTACATGATGTGGAGAGAGCCGAACAGGAGGAAGTGGAG
TGTGAGGAGAACGATGGGAGGAGGGCAGAGCGATGGCACCCACC
CGGTGGGAACCCCCCCCAGATAGCCCCGGAAATACACTGAACAGCTCACC
ACCCCTGGCTCCGGGAGTTGATCCGAACATCGACGTCTATGGAATT
AAGTGCCATGAAAACTCGCCTAGGAAGGAGGTGTACTTCATGGCAATTA
TTGACATCCTACTCATTATGATGCAAAAAGAAAGCTGCCATGCTGC
AAAAACTGTTAACATGGCGCTGGCGCGGAGATCTCCACCGTGAACCCA
GAACAGTATTCAAAGCGTTTGACTTTATTGGCACATCTGACGt
aagcggccgc